

CAN Module Upgrade

This guide explains the installation of the Barrett Technology CAN Module Upgrade unit.

Step 1: Turn OFF power to the WAM.


Step 2: Unlatch the four clips at the sides of the WAM base to remove the wrap-around base cover.

Step 3: Using a 4mm hex wrench, remove the flat socket-head cap screw at the upper left corner of the WAM back plate, and allow the back plate to flip down.




Step 4: Using a Phillips screwdriver, remove the two screws (circled in RED, above) holding the safety board cover in place.

Step 5: Identify the open CAN socket on the Safety Board. The CAN socket is circled in RED below, seen viewed from above and behind the WAM.

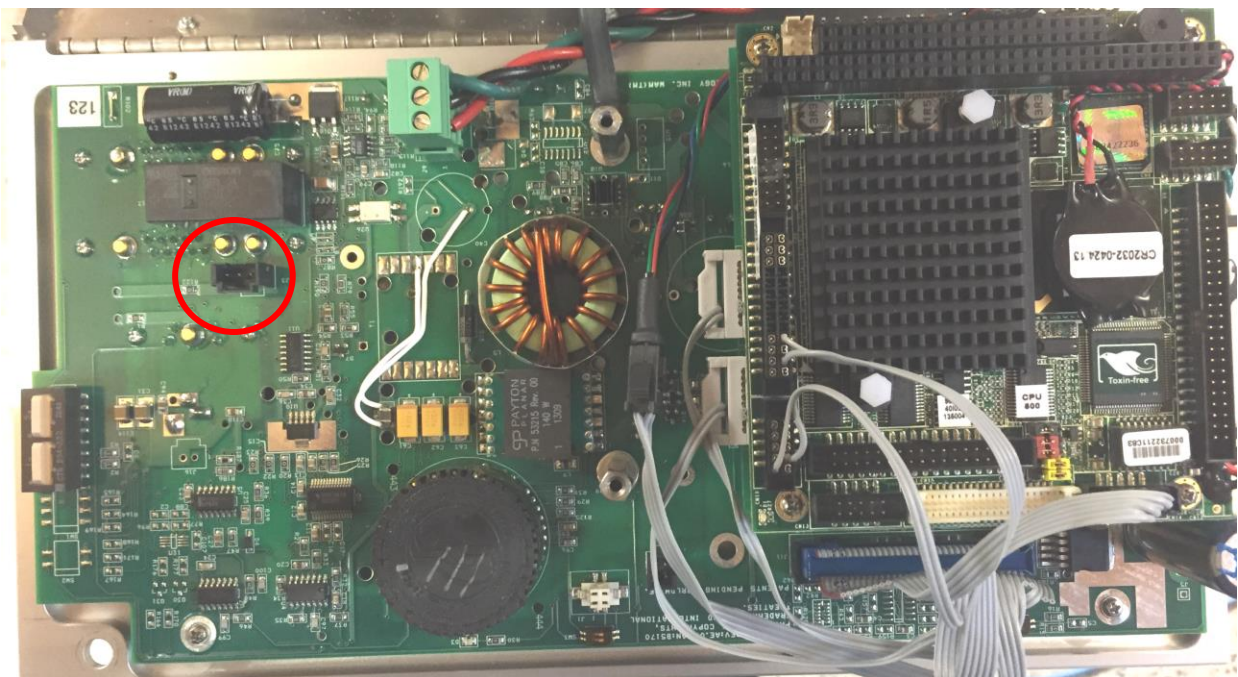

Step 6: Install the CAN Module Upgrade into the socket as shown below. Ensure that the locking tab on the Module Upgrade engages with the mating feature on the socket.

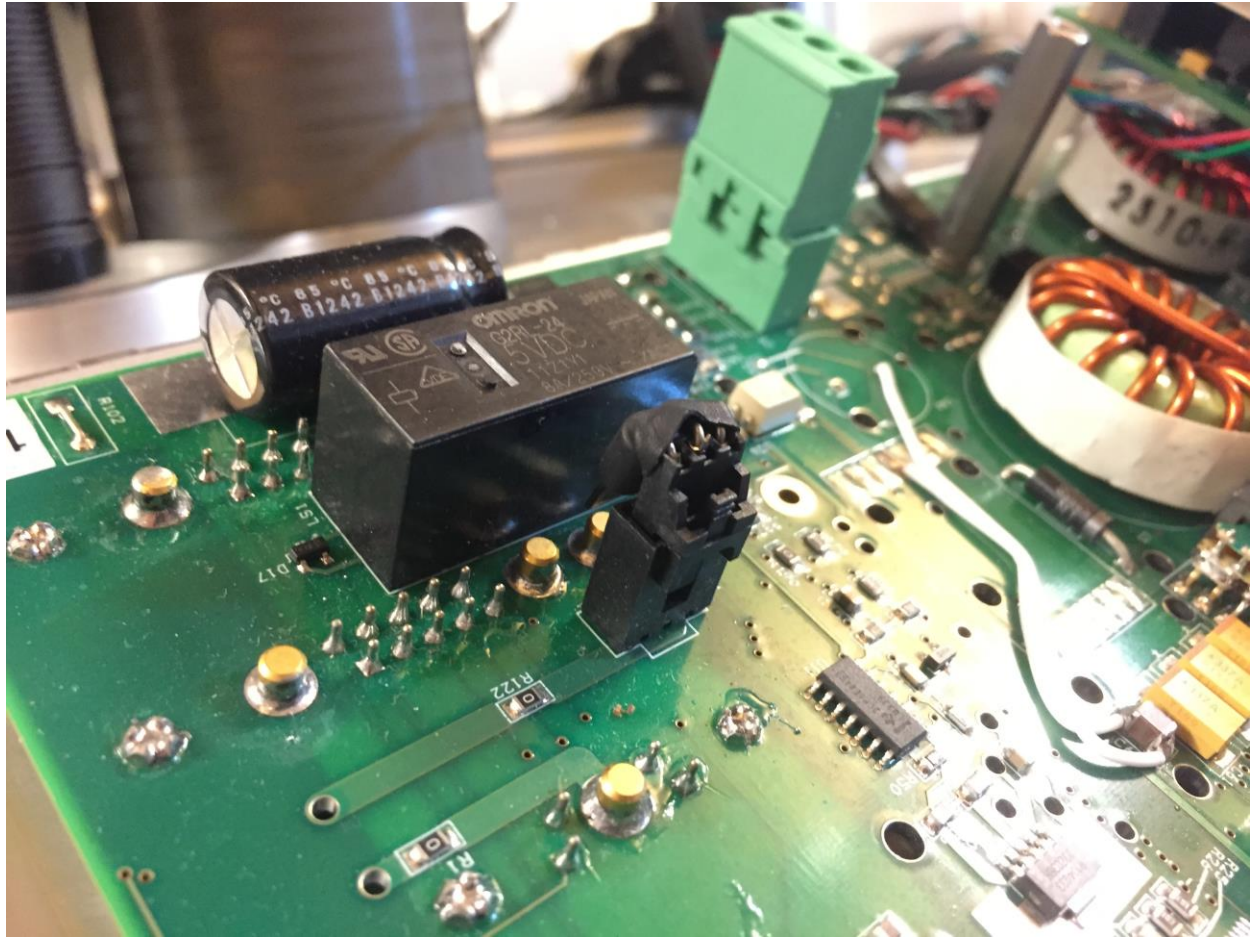

Step 7: Reinstall the safety board cover and close the WAM by reversing Steps 4, 3, and 2.

If you experience any difficulties during installation, please contact Barrett Technical Support at support@barrett.com.